

Pharmacy  
fridges  
2017/2018



**LIEBHERR**

Quality, Design and Innovation



# Pharmacy fridges compliant with DIN 58345

Fridges for use in pharmacies, hospitals and surgeries must meet very stringent requirements for the storage of pharmaceutical products. Liebherr pharmacy fridges compliant with DIN 58345 are therefore equipped with many functions and features to ensure optimum protection of high-quality preparations and sensitive pharmaceutical products. The flush-mounted electronic controller enables accurate temperature control and creates optimum storage conditions in conjunction with highly effective insulation and the forced-air cooling system. Integrated visual and audible alarm systems warn the user of undesired temperature deviations. Liebherr fridges offer the safety and consistency that are essential for the storage of pharmaceutical products: 24 hours a day, 365 days a year.

## Safety compliant with DIN 58345

All Liebherr medical storage appliances comply with the DIN standard "Fridges for drugs DIN 58345". This guarantees that the appliances are particularly suitable for medicine storage and that they offer the highest standard for conscientious users.

DIN 58345 includes the following basic requirements:

- Maintenance of an operating temperature from +2°C to +8°C
- Can be used in ambient temperatures from +10°C to +35°C
- Noise emission less than 60 dB(A)
- Visual and audible warning installation for high and low temperature alarms
- Preparation of operating temperature recording
- Visual and audible power failure alarm for at least 12h
- Safety thermostat to avoid temperatures below +2°C
- Mechanical resistance of internal fittings at least 100 kg/m<sup>2</sup>
- Lockable door

## Climate class definitions

Climate class SN: Ambient temperatures from +10°C to +32°C

Climate class N: Ambient temperatures from +16°C to +32°C

Climate class ST: Ambient temperatures from +16°C to +38°C

Climate class T: Ambient temperatures from +16°C to +43°C

# Good reasons to choose Liebherr



## Highest performance

Liebherr medical storage fridges offer constant refrigeration performance and maximum temperature consistency. The use of ultra-modern components, efficient and environmentally friendly refrigerants as well as precision controls ensures that pharmaceutical products and medicines are stored correctly. Optional monitoring software continuously logs the temperature profile – and if required, warning systems alert the user when defined temperature limits are exceeded.

## Efficiency

Highly efficient with low energy consumption: the precision electronic controller, the forced-air refrigeration system and highly effective insulation guarantee low operating costs – and are eco-friendly as well. The high quality of the appliances guarantees long life and reliable operation and provides for economic and ecological sustainability of medicine storage.

## Sturdy quality

Liebherr appliances are specially constructed for intensive professional use and feature a very sturdy design – with highest-quality materials and perfect workmanship right down to the smallest detail. The quality of the appliances is ensured by elaborate tests. Safety is given top priority in every pharmacy model. Many features ensure that pharmaceutical products are always protected – for example the sturdy wire shelves have a loading capacity of up to 45 kg.

## Easy cleaning

Hygiene plays a special role in medicine storage. That is why Liebherr appliances are equipped with moulded inner liners with large radii that are easy to clean. The shelves can easily be removed and cleaned. And the flush-mounted electronic controller with a membrane keypad is extremely resistant to soiling. 30 mm high roller sets can be retrofitted to the table-height MKUv models and adjustable legs can be retrofitted to the MKv models to allow easy cleaning underneath.

## Safety

Extremely safe and reliable: because stringent requirements with regard to temperature consistency must be met for the storage of pharmaceutical products, Liebherr appliances are equipped with many useful functions. A visual and audible alarm alerts the user to temperature deviations – and if the door is open for longer than 1 minute. The appliances are equipped with a volt-free contact for alarm forwarding to an external remote warning system. In addition, they have an RS 485 serial interface which enables as many as 20 appliances to be networked with a central documentation and alarm system.

## Easy servicing

Durability and reliability are given top priority in the material selection and development of Liebherr appliances. Their quality and functionality are ensured by elaborate tests. The LED ceiling lighting with a separate switch on glass door appliances provides for optimum illumination and allows fast access to the pharmaceutical products. To prevent unnecessary loss of cold air, the door is self-closing. The inner liners are designed with large corner radii for easy cleaning. Variable wire shelves enable easy and individual organisation of the interior.

# The advantages at a glance



Volt-free contact: For forwarding alarms to an **external remote warning system**.



**Visual and audible alarms** alert the user if the internal temperature deviates beyond +2°C and +8°C.

The **moulded inner liner** is exceptionally easy to clean and enables **perfect hygiene**.

The **evaporator cover** prevents contact between the **stored pharmaceutical products** and the cold evaporator surface.



The narrow-mesh **plastic-coated wire shelves** are easily adjustable with many shelf positions for maximum flexibility and use of space. The shelves are removable at 90° door opening.

Liebherr pharmacy fridges comply with **DIN 58345** for the storage of pharmaceutical products, thus ensuring the highest standards for protecting temperature-critical pharmacy products.



Access port (diameter 10 mm) allows **integration of an external temperature probe** (e.g. PT 100 or similar) in case independent temperature monitoring is required.



The precise **electronic controller** with digital display enables temperatures to be set accurately to 1/10°C. The large display allows improved temperature visibility at a glance.



The **LED ceiling light** with separate switch on glass door models MKv 3913 and MKUv 1613 allows good visibility of the fridge contents.



The flush-mounted controller has an **integrated data memory**. It logs e.g. the maximum and minimum internal temperatures that occur once the set temperature is reached. The min/max temperature memory function stores the values for up to 41 days.



Temperature is set at the factory to +5°C. The pharmacy fridges are equipped with an additional **safety thermostat** to prevent temperatures dropping below +2°C in case of a malfunction. In view of the potentially high value of the contents of a pharmacy fridge, this important feature gives added protection.

**Self-closing doors** for convenient use and improved temperature stability. The integrated lock is very sturdy and protects the stored products against unauthorised access.



In case of power failure, the electronic controller is immediately supplied by an **integrated 12 V battery** and indicates the failure optically and audibly for at least 12 hours.

MKv 3913

# Quality right down to the smallest detail

Liebherr medical storage fridges comply with DIN 58345. A visual and audible alarm alerts the user to temperature deviations or a power failure. The safety thermostat prevents temperatures below +2°C. All appliances are equipped with a volt-free contact

for alarm forwarding to an external remote warning system. The inner liner with the individually height-adjustable, sturdy wire shelves is easy to clean and enables perfect hygiene. This ensures safe storage of your pharmaceutical products.



#### Electronic controls.

The precision electronic controller features a 1/10°C temperature display. Symbols indicate the operating status of the appliance. To ensure hygiene, the electronic controller is flush-mounted and features a membrane keyboard. The temperature is set to +5°C at the factory. The pharmacy fridges are equipped with a safety thermostat to prevent temperatures below +2°C in case of a malfunction. The large display allows improved temperature visibility at a glance.



#### Integrated alarm systems.

Visual and audible alarm systems alert the user to undesired temperature deviations or when the door is open. All alarm parameters are pre-set in accordance with DIN 58345 requirements. The door-open alarm delay can be adjusted between 1 and 5 minutes. Additional safety is provided by a visual power failure alarm and sensor defect alarm.



#### Integrated data memory.

The flush-mounted electronic controller has an integrated data memory which logs the maximum and minimum internal temperatures as well as the last three temperature alarms and power failures, together with the time, date and duration. These data can be read out on the display.



#### External temperature sensor.

The pharmacy fridges have an access port (diameter 10 mm) on the rear which allows a temperature sensor to be positioned in the interior.

#### External temperature and alarm documentation.

The pharmacy fridges are equipped with a volt-free contact for alarm forwarding to an external remote warning system, as well as with a serial interface RS 485 which enables central documentation of temperature profiles and alarm events (optionally available).

#### Safety compliant with DIN 58345.

Liebherr medical storage fridges comply with the standard "DIN 58345". Therefore they are able to store pharmaceutical products at an operating temperature from +2°C to +8°C. Visual and audible alarm systems alert the user to undesired temperature deviations and power failures. A safety thermostat prevents temperatures below +2°C. In addition the appliances can be used in ambient temperatures from +10°C to +35°C – and can be locked to prevent unauthorised access.



#### Hygienic interior for versatile use.

Added hygiene: the seamless inner liner is exceptionally easy to clean. Moulded trayslides prevent the grid shelves from tipping and allow flexible height adjustment in increments of 32 mm. The evaporator cover prevents pharmaceutical products from freezing onto the rear wall.



#### Close-mesh wire shelves.

The plastic-coated shelves have an extremely high loading capacity of up to 45 kg. They are easily adjustable in height and can be removed at a door opening angle of 90°. Their close-mesh design ensures secure storage even for small items.



#### LED ceiling lighting.

The appliances with glass door (MKv 3913, MKUv 1613) have a LED ceiling light with separate switch. It ensures uniform illumination of the interior and enables fast, direct access to the stored pharmaceutical products



**Pharmacy fridges**  
compliant with DIN 58345

<b>Total gross / net capacity</b>	
<b>Exterior dimensions in mm (w/d/h)</b>	
<b>Interior dimensions in mm (w/d/h)</b>	
Energy consumption in 365 days <sup>1</sup>	
Climate rating	
Refrigerant	
Sound power level	
Voltage/Connection rating	
Cooling system / Defrost	
Temperature range	
External cabinet finish	
Door / lid material	
Internal liner material	
Type of control	
Temperature display	
Power failure alarm	
Fault: warning signal	
Interface / Volt-free alarm contact	
Interior light	
Adjustable shelves	
Usable shelf area in mm (w/d)	
Shelf material	
Shelf loading	
Handle	
Lock	
Self-closing door	
Door hinges	
Gross / net weight	
<b>Accessories</b>	
Shelf, plastic-coated	
Stacking kit, white	
Roller bars	
Adjustable legs	
RS 485/RS 232 adapter, data log software	
Additional lock barrels (up to 10)	

	<b>MKv 3913</b> MediLine	<b>MKv 3910</b> MediLine	<b>MKUv 1613</b> MediLine	<b>MKUv 1610</b> MediLine
	LED		LED	
Total gross / net capacity	360/281 litre	360/300 litre	141 / 115 litre	141 / 116 litre
Exterior dimensions in mm (w/d/h)	600/615/1840	600/615/1840	597/615/825	597/615/825
Interior dimensions in mm (w/d/h)	440/435/1635	440/435/1635	440/435/670	440/435/670
Energy consumption in 365 days <sup>1</sup>	480 kWh	309 kWh	369 kWh	273 kWh
Climate rating	SN	SN	SN	SN
Refrigerant	R 600a	R 600a	R 600a	R 600a
Sound power level	48 dB(A)	48 dB(A)	47 dB(A)	47 dB(A)
Voltage/Connection rating	220–240V~/1.5 A	220–240V~/1.5 A	220–240V~/1.0 A	220–240V~/1.0 A
Cooling system / Defrost	forced-air / automatic	forced-air / automatic	forced-air / automatic	forced-air / automatic
Temperature range	+5°C	+5°C	+5°C	+5°C
External cabinet finish	steel / white	steel / white	steel / white	steel / white
Door / lid material	glass door	steel	glass door	steel
Internal liner material	commercial grade moulded polystyrol	commercial grade moulded polystyrol	commercial grade moulded polystyrol	commercial grade moulded polystyrol
Type of control	electronic controller	electronic controller	electronic controller	electronic controller
Temperature display	external digital	external digital	external digital	external digital
Power failure alarm	immediately upon power failure for at least 12hrs	immediately upon power failure for at least 12hrs	immediately upon power failure for at least 12hrs	immediately upon power failure for at least 12hrs
Fault: warning signal	visual and audible	visual and audible	visual and audible	visual and audible
Interface / Volt-free alarm contact	RS 485 / yes	RS 485 / yes	RS 485 / yes	RS 485 / yes
Interior light	LED lighting, with separate switch		LED lighting, with separate switch	
Adjustable shelves	5	5	3	3
Usable shelf area in mm (w/d)	440 / 420	440 / 420	440 / 420	440 / 420
Shelf material	wire shelves plastic-coated	wire shelves plastic-coated	wire shelves plastic-coated	wire shelves plastic-coated
Shelf loading	45 kg	45 kg	45 kg	45 kg
Handle	ergonomic slimline handle	ergonomic slimline handle	ergonomic slimline handle	ergonomic slimline handle
Lock	fitted	fitted	fitted	fitted
Self-closing door	yes	yes	yes	yes
Door hinges	right, reversible on site	right, reversible on site	right, reversible on site	right, reversible on site
Gross / net weight	91 / 85 kg	74 / 68 kg	47 / 45 kg	41 / 39 kg
<b>Accessories</b>				
Shelf, plastic-coated	7112313	7112313	7112313	7112313
Stacking kit, white			9876687	9876687
Roller bars			9590521	9590521
Adjustable legs	9590231	9590231		
RS 485/RS 232 adapter, data log software	9590387	9590387	9590387	9590387
Additional lock barrels (up to 10)	on request	on request	on request	on request

<sup>1</sup> Measured at +25°C ambient temperature and +5°C set temperature

Shelves plastic-coated

Additional plastic-coated shelves can be retrofitted as required to provide extra storage options. They have an extremely high loading capacity of up to 45 kg.



Stacking kit

A stacking kit is available as an accessory for combining the MKUv 1613 and MKUv 1610 models as well as other table-height appliances in the Liebherr research and laboratory range. This allows stacking a glass door appliance and solid door appliance or a fridge and freezer on a small footprint.



Roller bars and adjustable legs

30 mm high roller bars can be retrofitted to the table-height MKUv models and adjustable legs can be retrofitted to the MKv models to allow easy cleaning underneath. The legs are adjustable between 115 mm and 170 mm in height.



Interface converter with monitoring software

A special interface converter including LTM monitoring software is available for central documentation of the temperature profile data and alarm events of multiple appliances via the RS 485 serial interface. Capability is provided for interconnecting as many as 20 appliances and centrally documenting their parameters. Alarm forwarding or regular status reports to up to three email addresses can be configured as an additional feature. Commercially available converters can also be used to connect the connected appliances to a PC via WLAN or LAN. Software requirements: PC with Windows® operating system.



Additional lock barrels

Up to 10 additional lock barrels to protect the stored products against unauthorised access are available as accessories for the pharmacy fridges. In case of multiple appliances, different locks can be installed to ensure that the respective appliance can only be accessed by the staff members responsible for it.



# Liebherr commercial appliances can be found at specialist retailers where providing first-class service and expert advice is their top priority.

You can find an overview of all Liebherr appliances in our main catalogues. Available in stores or on [home.liebherr.com](http://home.liebherr.com).



## Smart communication on all fronts



Discover the world of Liebherr-Appliances with the latest interesting news, stories, valuable tips and tricks on food storage, recipes and lots more besides!



### WineGuide App

A brief overview of Bordeaux wines and their producers, plus storage information.



[apps.home.liebherr.com](http://apps.home.liebherr.com)

Find out which app is available for which operating system (Apple, Android etc.), and which device type.



The Liebherr-Hausgeräte YouTube channel provides interesting and helpful films showing you the functions of Liebherr fridges and freezers.



News, product launches and special promotions can be found on our Facebook page, the Liebherr blog, on Instagram and Pinterest.



[socialmedia.home.liebherr.com](http://socialmedia.home.liebherr.com)

Learn all about our current social media channels.



Subject to modification. Current data see [home.liebherr.com](http://home.liebherr.com). Printed in Germany by Typodruck 7944986-02/2.5/10.2017